

细节见证品质
Details define quality.



品质成就未来
Quality makes future.

振动筛分机
Vibrating Screen Powder Machine

使用说明书

OPERATING MANUAL

<使用前请仔细阅读>
Please read this manual carefully before use.

兰博涂装科技有限公司
LANBO COATING TECHNOLOGY CO.,LTD.

WWW.LVOBO.COM

400 605 9596




兰博涂装科技(泉州)有限公司
LANBO COATING TECHNOLOGY CO.,LTD

LVOBO[®]

科技 | 创新 | 品质 | 服务

» BO-200SY (加上下振幅款) 实验筛

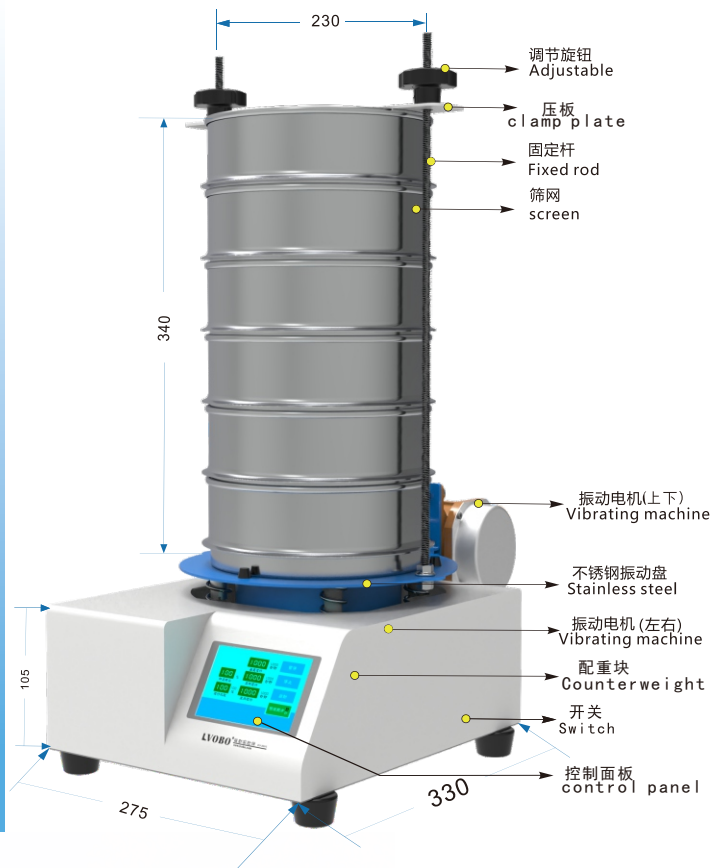

操作说明



1. 按下背部电源总开关等待屏幕亮起
2. 该设备具有三种运行方式(垂直 正转 反转)，若三种方式均有设定时间(≤1000S)，则物料先正转运行，运行时间结束后垂直运行，再反转后垂直(正转——垂直——反转——垂直)，垂直运行作用于物料卡在筛网时的上下震动
3. 可以根据所需调整转数输出
4. 运行次数根据所需设定次数(最高100次)，每种运行停止为一次
5. 设定数值后，点击启动设备开始运行，若想中断点击暂停，从头运行点击停止
6. 可根据物料种类设定不同运行方式保存到预设配方

技术参数

PARAMETER

| | | | |
|--------------------------------------|--|-------------------------------------|-------------------------------------|
| 产品型号 Product model | BO-200SY(加上下震动) BO-200SY(常规款) | 振动频率 Frequency of vibration | 1400(n/MIN) |
| 输入电压(可选) Input voltage (optional) | AC220V 50Hz 单相 | 激振功率 Vibration rate | 1(KN) |
| 有效筛面积 Effective screening area | Φ200*H100(MM) | 电机功率 Power of motor | 200W |
| 筛网层数 Screen layers | 1-6层(每层为45mm) 1-6layer (Each layer is 45mm) | 不锈钢参数 Stainless steel parameters | SUS304不锈钢 SUS304 Stainless steel |
| 筛网直径 Screen diameter | Φ200(MM) | 筛网目数 Screen number | 2-2000目(可选) 2-2000eyes(Optional) |

注意事项 Cautions

- 一、使用设备前，请检查设备是否有损伤。
Please check the device for any damage prior to use.
- 二、连接电源前，请检查设备开关是否处于关闭状态。
Please make sure the start switch is turned off before connecting to the power supply.
- 三、将密封条装置在筛斗边缘上，正确装上筛网，用弹簧挂钩加以固定。
Please install the sealing strip on the edge of the screen bucket and fix with the spring hook.
- 四、将所需筛分的产品倒入筛网内，机台经过电机激振，产品往筛斗出料口处流出即可。
Please pour the required products into the screen, after working, the end products will flow out from the outlet port.
- 五、注意清洁筛斗，增加筛斗使用寿命。
Please clean the screen bucket after use to increase the service life of the device.

操作说明 operating instructions

- 一、本说明书提及的安全事项必须阅读和理解，然后才能使用设备。
The mentioned cautions must be read and understood before using the device.
- 二、设备不得暴露在雨天和潮湿的环境中，必须处于常温干燥环境中使用；且远离可燃性气体、液体。
The device shall not be exposed to rainy and wet environment, it must be used in normal temperature dry environment, and stay away from combustible gases and liquids.
- 三、使用设备前，确保安全和正常工作。
Please ensure the safe and normal operation before using the device.
- 四、每次使用时必须检查输入电源电压是否合格。并接入符合参数标准的插头。
It is necessary to check whether the input power voltage is qualified during each use. And connect the plug that meets the parameter standards.
- 五、设备必须对大地连接
The equipment must be connected to the earth.

